

KAMIYA BIOMEDICAL COMPANY

Human Beta 2-Microglobulin ELISA

**For the quantitative determination of Beta 2-Microglobulin
in human biological fluids**

Cat. No. KT-500

For Research Use Only. Not for use in diagnostic procedures in the U.S.

PRODUCT INFORMATION**Human Beta 2-Microglobulin ELISA**
Cat. No. KT-500**INTENDED USE**

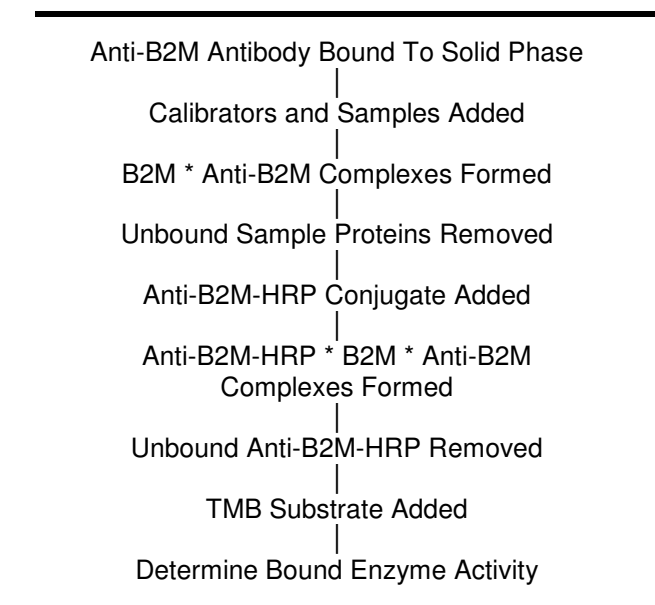
The Human Beta 2-Microglobulin ELISA is a highly sensitive two-site enzyme-linked immunoassay (ELISA) for the quantitative determination of Beta 2-Microglobulin in human biological fluids. For research use only.

INTRODUCTION

Beta 2-Microglobulin (B2M) is an 11 kDA protein. It forms the subunit of the MHC class I molecule and associates with the outer membrane of many cells including lymphocytes. It is present in low levels in serum and urine of normal people, but at a higher concentration in people with renal diseases, kidney transplants and various other inflammatory and infectious conditions.

PRINCIPLE

The principle of the double antibody sandwich ELISA is represented in Figure 1. In this assay the Beta 2-Microglobulin present in samples reacts with the anti-Beta 2-Microglobulin antibodies which have been adsorbed to the surface of polystyrene microtiter wells. After the removal of unbound proteins by washing, anti-B2M antibodies conjugated with horseradish peroxidase (HRP), are added. These enzyme-labeled antibodies form complexes with the previously bound B2M. Following another washing step, the enzyme bound to the immunosorbent is assayed by the addition of a chromogenic substrate, 3,3',5,5'-tetramethylbenzidine (TMB). The quantity of bound enzyme varies directly with the concentration of B2M in the sample tested; thus, the absorbance, at 450 nm, is a measure of the concentration of B2M in the test sample. The quantity of B2M in the test sample can be interpolated from the calibration curve constructed from the calibrators, and corrected for sample dilution.

Figure 1.**COMPONENTS**

1. Diluent Concentrate
One bottle containing 50 mL of a 5X concentrated diluent running buffer.
2. Wash Solution Concentrate
One bottle containing 50 mL of a 20X concentrated wash solution.

3. Enzyme-Antibody Conjugate Concentrate
One vial containing 150 μ L of a 100X concentrated affinity-purified anti-human Beta 2-Microglobulin antibody conjugated with HRP in stabilizing buffer.
4. TMB Substrate Solution
One vial containing 12 mL of TMB and hydrogen peroxide in citric acid buffer at pH 3.3.
5. Stop Solution
One vial containing 12 mL of 0.3 M sulfuric acid. WARNING: Avoid contact with skin.
6. Microtiter Plate
Twelve removable eight-well strips in well holder frame. Wells are coated with affinity-purified anti-human B2M.
7. Human Beta 2-Microglobulin Calibrator
One vial containing a lyophilized Human Beta 2-Microglobulin Calibrator.
8. Positive Control
One vial containing 50 μ L of serum with 0.1% sodium azide. See the Control Certificate for the concentration.

MATERIALS REQUIRED BUT NOT PROVIDED

- Precision pipettes (2 μ L to 200 μ L) for making and dispensing dilutions
- Test tubes
- Microplate washer/aspirator
- Distilled or de-ionized H₂O
- Microplate reader
- Assorted glassware for the preparation of reagents and buffer solutions
- Timer

PRECAUTIONS

1. Read the instructions carefully before beginning the assay.
2. This kit is for research use only.
3. Great care has been taken to ensure the quality and reliability of this product. However, it is possible that in certain cases, unusual results may be obtained due to high levels of interfering factors.
4. Preservatives
 - Positive Control contains 0.1% sodium azide.
5. No additives or preservatives are necessary to maintain the integrity of the specimen. Avoid azide contamination.
6. Azide and thimerosal at concentrations higher than 0.1% inhibit the enzyme reaction.
7. Other precautions:
 - Do not interchange kit components from different lots.
 - Do not use kit components beyond the expiration date.
 - Protect reagents from direct sunlight.
 - Do not pipette by mouth.
 - Do not eat, drink, smoke or apply cosmetics where reagents are used.
 - Avoid all contact with the reagents by using gloves.
 - Stop solution contains diluted sulfuric acid. Irritation to eyes and skin is possible. Flush with water after contact.

REAGENT PREPARATION

1. Diluent Concentrate
The Diluent solution supplied is a 5X concentrate and must be diluted 1:5 with distilled or de-ionized water.
2. Wash Solution Concentrate
The Wash Solution supplied is a 20X concentrate and must be diluted 1:20 with distilled or de-ionized water. Crystal formation in the concentrate is not uncommon when storage temperatures are low. Warming of the concentrate to 30-35°C before dilution can dissolve crystals.

3. Enzyme-Antibody Conjugate Concentrate
Calculate the required amount of working conjugate solution for each microtiter plate test strip by adding 10 μL Enzyme-Antibody Conjugate to 990 μL of 1X Diluent for each test strip to be used for testing. Mix uniformly, but gently. Avoid foaming.
4. TMB Substrate Solution
Ready to use as supplied.
5. Stop Solution
Ready to use as supplied.
6. Microtiter Plate
Ready to use as supplied. Unseal microtiter pouch and remove plate from pouch. Remove all strips and wells that will not be used in the assay and place back in pouch and re-seal along with desiccant.
7. Human Beta 2-Microglobulin Calibrator
Add 1.0 mL of distilled or de-ionized water to the lyophilized Human B2M Calibrator and mix gently until dissolved. The calibrator is now at a concentration of 710 ng/mL (the reconstituted calibrator should be aliquoted and frozen if future use is intended). Prepare the Human B2M Calibrators immediately prior to use according to the table below. Mix well between each step. Avoid foaming.

Calibrator	Concentration (ng/mL)	Calibrator Volume added to 1X Diluent	Volume of 1X Diluent
6	100	100 μL Human B2M Calibrator	610 μL
5	50	300 μL Calibrator 6	300 μL
4	25	300 μL Calibrator 5	300 μL
3	12.5	300 μL Calibrator 4	300 μL
2	6.25	300 μL Calibrator 3	300 μL
1	3.13	300 μL Calibrator 2	300 μL
0	0		600 μL

8. Positive Control
The concentration and recommended dilution are provided on the Control Certificate. Before use, briefly centrifuge the Positive Control to allow all of the liquid to collect in the bottom of the vial.

STORAGE AND STABILITY

1. Complete Kit
The expiration date for the kit is stated on the outer label. The recommended storage temperature is 4°C. **Note: See long term storage recommendations below for the Human Beta 2-Microglobulin Calibrator and Positive Control.**
2. Diluent
The 5X Diluent Concentrate is stable until the expiration date. The 1X working solution is stable for at least one week from the date of preparation. Both solutions should be stored at 4°C.
3. Wash Solution
The 20X Wash Solution Concentrate is stable until the expiration date. The 1X working solution is stable for at least one week from the date of preparation. Both solutions can be stored at room temperature (RT, 16-25°C) or at 4°C.
4. Enzyme-Antibody Conjugate
Undiluted anti-B2M-HRP conjugate should be stored at 4°C and diluted immediately prior to use. The working conjugate solution is stable for up to 1 hour when stored in the dark.
5. TMB Substrate Solution
The TMB Substrate Solution should be stored at 4°C and is stable until the expiration date.

6. Stop Solution
The Stop Solution should be stored at 4°C and is stable until the expiration date.
7. Microtiter Plate
Anti-human B2M coated wells are stable until the expiration date, and should be stored at 4°C in the sealed foil pouch with desiccant pack.
8. Human Beta 2-Microglobulin Calibrator
The lyophilized Human Beta 2-Microglobulin Calibrator should be stored at 4°C or frozen until reconstituted. The reconstituted calibrator should be aliquoted and stored frozen. Avoid multiple freeze/thaw cycles. The working calibrator solutions should be prepared immediately prior to use and are stable for up to 1 hour.
9. Positive Control
For storage longer than 7 days keep frozen until the expiration date. Storage less than 7 days can be at 4°C. Avoid multiple freeze/thaw cycles.

INDICATIONS OF INSTABILITY

If the test is performing correctly, the results observed with the calibrator solutions should be within 20% of the expected values.

SPECIMEN COLLECTION AND HANDLING

Blood should be collected by venipuncture and the serum separated from the cells, after clot formation, by centrifugation. For plasma samples, blood should be collected into a container with an anticoagulant and then centrifuged. Care should be taken to minimize hemolysis, excessive hemolysis can impact your results. Assay immediately or aliquot and store samples at -20°C. Avoid repeated freezing/thawing.

For any sample that might contain pathogens, care must be taken to prevent contact with open wounds. No additives or preservatives are necessary to maintain the integrity of the specimen. Avoid azide contamination.

ASSAY PROTOCOL

Dilution of Samples

The assay for quantification of B2M in samples requires that each test sample be diluted before use. For a single step determination a dilution of 1:200 is appropriate for most serum/plasma samples. For absolute quantification of samples that yield results outside the range of the calibration curve, a lesser or greater dilution might be required. If unsure of sample level, a serial dilution with one or two representative samples before running the entire plate is highly recommended.

To prepare a 1:200 dilution of sample, transfer 4 µL of sample to 796 µL of 1X Diluent. This gives you a 1:200 dilution of your sample. Mix thoroughly.

Procedure

Bring all reagents to RT before use.

1. Pipette 100 µL of
 - Calibrator 0 (0.0 ng/mL) in duplicate
 - Calibrator 1 (3.13 ng/mL) in duplicate
 - Calibrator 2 (6.25 ng/mL) in duplicate
 - Calibrator 3 (12.5 ng/mL) in duplicate
 - Calibrator 4 (25 ng/mL) in duplicate
 - Calibrator 5 (50 ng/mL) in duplicate
 - Calibrator 6 (100 ng/mL) in duplicate
2. Pipette 100 µL of Positive Control (in duplicate) into pre-designated wells.
3. Pipette 100 µL of diluted sample (in duplicate) into pre-designated wells.

4. Incubate the Microtiter Plate at 22°C (RT) for sixty (60 ± 2) minutes. Keep plate covered and level during incubation.
5. Following incubation, aspirate the contents of the wells.
6. Completely fill each well with appropriately diluted Wash Solution and aspirate. Repeat three times, for a total of four washes. If washing manually: completely fill wells with diluted Wash Solution, invert the plate and pour/shake out the contents in a waste container. Follow this by sharply striking the wells on absorbent paper to remove residual Wash Solution. Repeat three times for a total of four washes.
7. Pipette 100 µL of appropriately diluted Enzyme-Antibody Conjugate to each well. Incubate at 22°C (RT) for sixty (60 ± 2) minutes. Keep plate covered in the dark and level during incubation.
8. Wash and blot the wells as described in Steps 5 and 6.
9. Pipette 100 µL of TMB Substrate Solution into each well.
10. Incubate in the dark at RT for precisely ten (10) minutes.
11. After ten minutes, add 100 µL of Stop Solution to each well.
12. Determine the absorbance at 450 nm of the contents of each well. Zero the plate reader to air.

The absorbance of the final reaction mixture can be measured up to 2 hours after the addition of the Stop Solution. However, good laboratory practice dictates that the measurement be made as soon as possible.

RESULTS

1. Subtract the average background value from the test values for each sample.
2. Using the results observed for the calibrators construct a calibration curve. The appropriate curve fit is that of a four-parameter logistics curve, although a second order polynomial (quadratic) or other curve fits may also be used.
3. Interpolate test sample values from calibration curve. Correct for sample dilution factor to arrive at Beta 2-Microglobulin concentration in original sample.

PERFORMANCE CHARACTERISTICS

In accord with good laboratory practice, the assays for specific Beta 2-Microglobulin require meticulous quality control. Each laboratory should use routine quality control procedures to establish inter- and intra-assay precision and performance characteristics.

LIMITATION OF THE PROCEDURE

1. Reliable and reproducible results will be obtained when the assay procedure is carried out with a complete understanding of the information contained in the package insert instructions and with adherence to good laboratory practice.
2. Factors that might affect the performance of the assay include proper instrument function, cleanliness of glassware, quality of distilled or de-ionized water, and accuracy of reagent and sample pipettings, washing technique, incubation time or temperature.

FOR RESEARCH USE ONLY

KAMIYA BIOMEDICAL COMPANY

12779 Gateway Drive, Seattle, WA 98168

Tel: (206) 575-8068 Fax: (206) 575-8094

Email: LifeScience@k-assay.com

www.k-assay.com